Ohaus Corporation
 29 Hanover Road
 Florham Park NJ
 07932-0900

Triple Beam Balance

Balanza de Triple Brazo

Instuction Manual

Please read this manual before you use your Ohaus Triple Beam Balance.

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Unpacking:

Carefully remove the balance and the separate poise (sliding weight) from the protective carton. You will find a slit rubber washer lodged underneath the platform, and one rubber washer located above the pointer. The washers are to be removed from the scale.

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Set-up:

After placing the balance on a smooth, flat surface, slide the separate poise up into the slot on the back of the center beam. Tilt poise over into place on the beam. With all poises in zero position, the pointer should be near zero. On applicable models, the tare poise (13) shall be at the extreme left of its bar.

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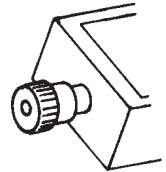
Zeroing:

For exact zero, adjust the knurled knob which is located at the left end of the beam. It is advisable to check the zero adjustment periodically.

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
Weighing:

Place the specimen on the center of the platform and proceed as follows:

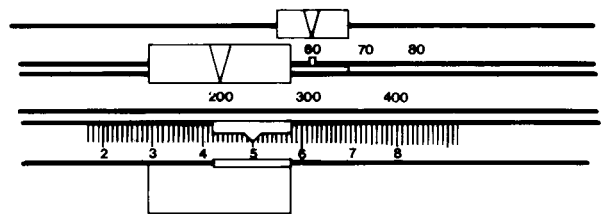1. Starting with the largest capacity beam (500 g), move the 500 g poise to the right to the first notch which causes the pointer to drop, then, move it back one notch, causing the pointer to rise.
2. Repeat procedure with the 100 g poise.
3. Slide the 10 g poise to the position which brings the pointer to rest at zero. The weight of the specimen is the sum of the values of all poise positions, read directly from the graduated beams.

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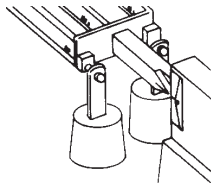
Attachment weights:

Total capacity is either 2610 grams or 5 pounds, 2 ounces when attachment weights are suspended from the pivots (14). Without the weights, the capacity is either 610 grams or 1 pound, 2 ounces.

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
Use of the tare:

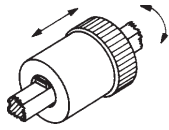Certain models are equipped with a patented tare poise (13). The poise will counterbalance empty containers no heavier than 225 grams or 8 ounces, by sliding it to the approximate balance, then rotating in either direction for precise positioning. Net weight of the contents of the container may then be read directly in the usual manner.

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Care and maintenance:

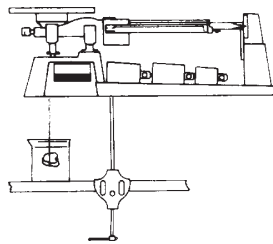
Keep balance clean at all times. In general, most foreign matter may be easily removed by an air syringe, but a piece of adhesive-backed tape pressed against the magnet faces will keep them free from dirt. Never apply lubricants to knives or bearings, nor allow foreign matter to accumulate.

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Specific gravity determination:

The Model 183 Clamp and Rod Support is an accessory for elevating the balance for suspending specimens in water. The rod is inserted into the 1/2 inch (12.7 mm) recess on the underside of the base. By means of a fine wire or thread, the specimen is attached to the check pin which connects the check rod assembly to the platform loop under the base.

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Formulas:

Sólidos - (denser than water): To determine specific gravity of solids denser than water, the specimen should be first weighed in air and then weighed immersed in water.

$$\frac{\text{Weight in air}}{\text{Weight in air minus weight in water}}$$

Sólidos - (less dense than water): To determine the specific gravity of solids less dense than water, it is necessary to attach a sinker and then make the following weighings:

- a. Weight of body in air.
- b. Combined weight of body in air plus sinker in water.
- c. Weight of both body and sinker in water.

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$$\frac{a}{b - c}$$

Features:

- A low cost, maintenance-free balance which provides all the convenience of a top loader, yet retains the ruggedness needed for a wide range of lab work.
- Tiered beam visibility and zero adjust combine to give high speed performance and reduce error.
- Magnetic damping speeds up weighing by causing the beam to come to rest quickly without affecting sensitivity or accuracy. It operates on the principle of a permanent magnetic field resisting the motion of a non-magnetic, aluminum damper vane attached to the beam. The pole faces of the damping magnets are positioned on both sides of the damper vane. Damping force is proportional to vane velocity and reduces to zero when the beam stops, thereby turning itself off. The system is permanent, self-regulating, maintenance-free, frictionless and effective at all loads.
- Tare beam and poise (13) (on specific models) make repetitive weighings and chemical formulations easy by eliminating container weight from calculations.
- Attachment weights store conveniently in the base when not in use.

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
Accessories

183-00 Specific Gravity Rod and Clamp
 703-00 Footed Polypropylene Scoop, 12" x 6" x 2-3/4" and Counterweight
 703-S0 Footed Stainless Steel Scoop, 12" x 6" x 2-3/4" and Counterweight
 706-00 Vinyl Dust Cover
 707-00 Metric Attachment Weight Set: two 1000g, one 500g
 807-03 Avoirdupoise Attachment Weight Set: two 2 lb, one 1 lb

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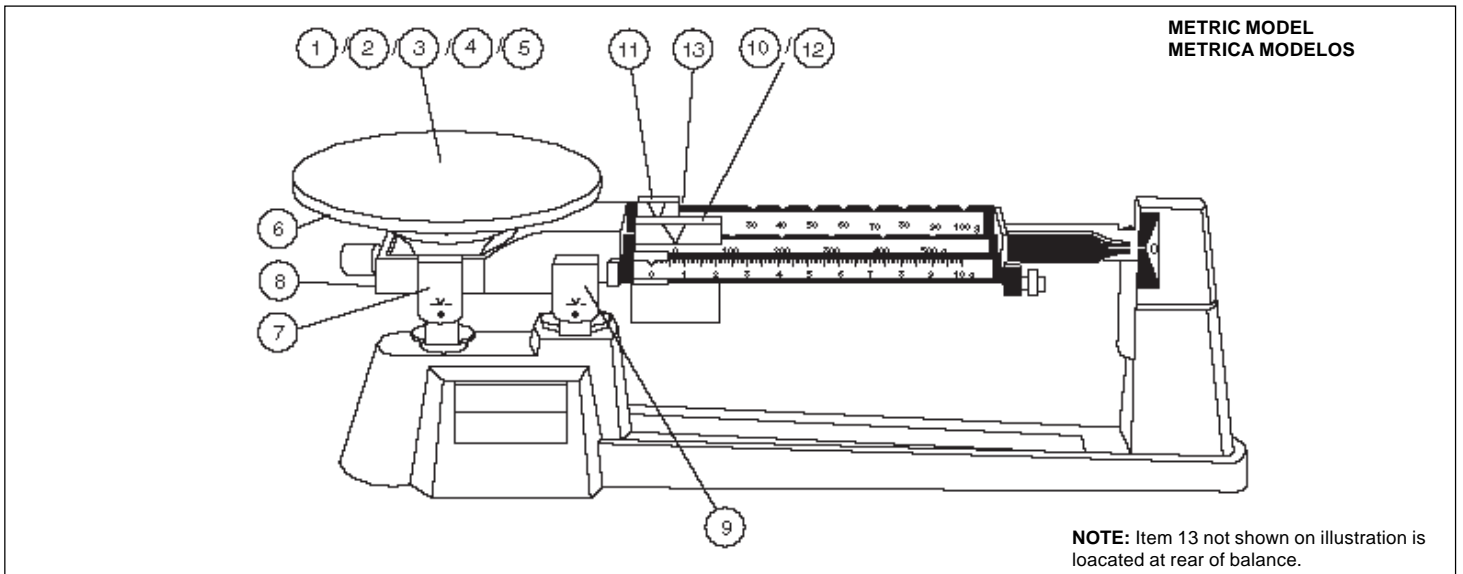Specifications

	700 Metric Series	800 Avoirdupois Series
Capacity		
w/attachment weights	2610g	5 lb. 2 oz.
w/o attachment weights	610g	1 lb, 2 oz.
Readability		
	0.1g	0.01 oz.
Calibrations		
Front Beam	10g x 0.1g	1oz. x 1/64 oz.
Center Beam	500g x 100g	16 oz. x 1 oz.
Rear Beam	100g x 10g	1oz. x 0.01 oz.

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**LIMITED WARRANTY**

Ohaus products are warranted against defects in, materials and workmanship from the date of delivery through the duration of the warranty period. During the warranty period Ohaus will repair, or, at its option, replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus.

This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product or as a result of service or modification by other than Ohaus. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages. As warranty legislation differs from state to state and country to country, please contact Ohaus or your local Ohaus dealer for further details.

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1600 Series Balance

Instuction Manual

Please read this manual before
unpacking, setting up and using
your balance.

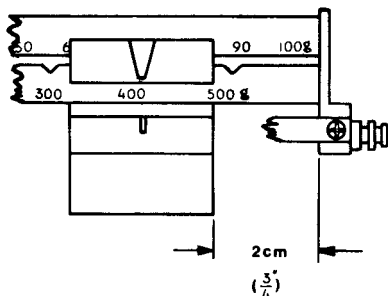
UNPACKING

Your new precision balance was protected during shipment by a specially designed carton. Carefully remove the balance, the attachment weights (included with some models) and the separately packaged Poise (sliding weight) from the carton and place the balance on a reasonably flat and level surface.


Remove the slit rubber washer, which is lodged underneath the platform support, and the plain rubber washer that is located above the pointer.

SET UP

Locate the bevel on the back of the center beam and slide the separate Poise up into it. Tilt the Poise over into place on the beam.

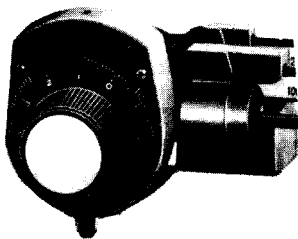Place both sliding Poises and the Dial at their zero positions.


Move the Tare Poise to the extreme left.

The Pointer should come to rest near the zero balance position.

ZEROING

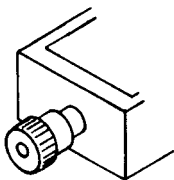To obtain exact zero balance, adjust the Balance Compensator (knurled knob) which is located at the left end of the beam.


It is recommended that the zero balance be checked periodically to insure accurate weight readings.

WEIGHING

When exact zero balance has been obtained, proceed as follows:

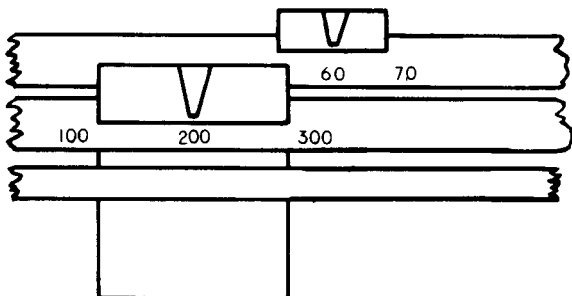1. Place the specimen in the center of the Platform.
2. Starting with the largest capacity beam (500 gram), move the 500 gram Poise toward the right to the first notch that causes the Pointer to drop, then move it back one notch, which causes the Pointer to rise.
3. Repeat the above procedure with the 100 gram Poise.


4. Turn the Dial until the Pointer comes to rest at zero.

The weight of the specimen is the sum of the values of the Poise positions, read directly from the Graduated Beams, and the position of the Dial.

OPTIONAL ATTACHMENT WEIGHTS

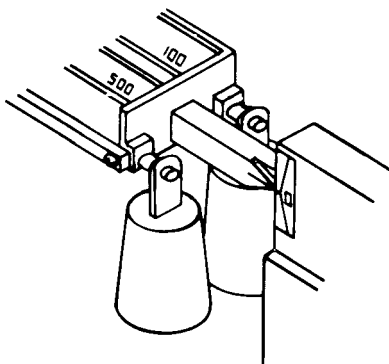The total capacity of the balance when Attachment Weights are suspended from the pivots is 2610 grams.


Without attachment weights, the capacity is 610 grams.

USE OF THE TARE

All models are equipped with a patented Tare Poise which will counterbalance empty containers up to and including 225 grams.

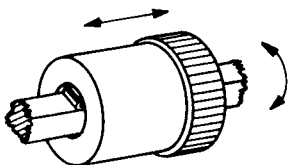To use the Tare Poise, slide it to a position of approximate balance, and then rotate the Poise to obtain exact zero balance.


The net weight of the contents of the container may then be determined in the usual manner.

FEATURES

- **CONVENIENT DIAL READING**

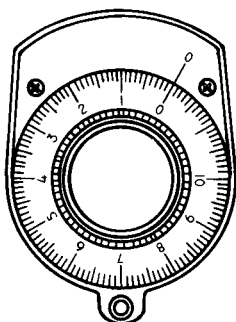Enables weighing to 0.1 gram without the need for an added poise.

- **TIERED BEAM VISIBILITY AND ZERO ADJUST**

Combination provides high speed efficiency and reduces errors.

- **MAGNETIC DAMPING**

Based on the principle of a permanent magnetic field resisting the motion of the non-magnetic aluminum damper vane. The damping force is proportional to the velocity of the damper vane and reduces to zero when the vane comes to rest.

Speeds up weighing by causing the beam to come to rest quickly without affecting sensitivity or accuracy. The system is permanent, self-regulating, frictionless and effective at all loads.

- **TARE BEAM AND POISE**

Simplifies repetitive weighings and chemical formulations by eliminating the container weight from calculations.

- **OPTIONAL ATTACHMENT WEIGHTS**

Store conveniently in the base when not in use.

CARE AND MAINTENANCE

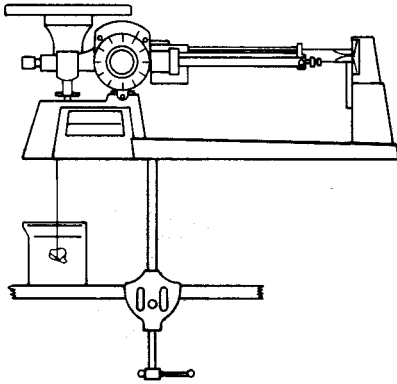
Keep the balance clean at all times. In general, most foreign matter may be easily removed by using an air syringe. A piece of adhesive-backed tape pressed against the Magnet faces will keep them free from dirt.

Never apply lubricants to the knives or bearings, nor allow foreign matter to accumulate.

SPECIFIC GRAVITY DETERMINATION

Ohaus Model 183 Clamp and Rod Support is an accessory which elevates the balance for suspending specimens in water.

The rod is inserted into the half inch (12.7 mm) diameter recess on the underside of the base. The specimen is attached to the check pin by means of a fine wire or thread. The check pin connects the check rod assembly to the platform loop under the base.

FORMULAS

- Solids denser than water.

First weigh the specimen in air and then weigh it immersed in water.

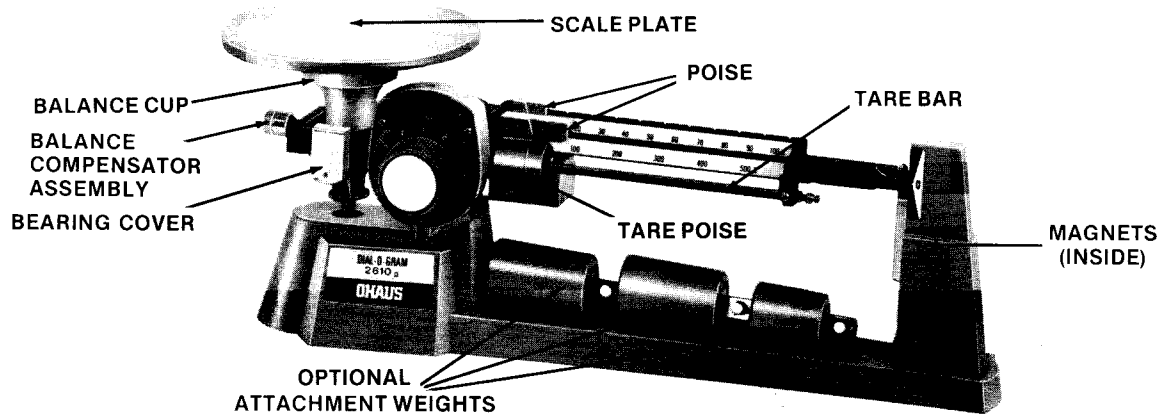
$$SG = \frac{\text{Weight in air}}{\text{Weight in air minus weight in water}}$$

- Solids less dense than water.

It is necessary to attach a sinker and then make the following weighings:

- a. Weight of specimen in air
- b. Combined weight of specimen in air plus sinker in water
- c. Weight of both specimen and sinker in water

$$SG = \frac{a}{b - c}$$

Model 1650 shown

REPLACEMENT PARTS LIST

Part No.	Description	Reqd	Models Used On
3045-00	Scale Plate, Stainless Steel 6" dia.	1	1650
1039-02	Scale Pan 6" dia. X 3/4"	1	1610
1064-30	Cross	1	
1077-03	Scoop, Stainless Steel 12" X 6" X 2 3/4" deep	1	1620-S
1101-20	Scoop, Polypropylene 12" X 6" X 2 3/4" deep	1	1620
3043-20	Specimen Pan	1	1630
3052-00	Cover	1	
2570-21	Cross	1	
1023-00	Balance Cup	1	All
1034-00	Friction Plate	3	All
1247-00	Friction Plate (with tab)	1	All
1008-05	Bearing	4	All
1035-03	Bearing Cover	4	All
3134-02	500g Poise Assembly	1	All
3550-04	Dial Housing Assembly	1	All
9701-01	Balance Compensator Assembly	1	All
9706-01	Tare Beam Assembly	1	All

ACCESSORIES

- 183-00 Specific Gravity Rod and Clamp
- 703-00 Footed Polypropylene Scoop,
12" X 6" X 2 3/4" and Counterweight
- 703-S0 Footed Stainless Steel Scoop,
12" X 6" X 2 3/4" and Counterweight
- 706-00 Vinyl Dust Cover
- 707-00 Metric Attachment Weight Set: two P/N 702-02,
1000g and one P/N 701-02, 500g
- 1078-03 Footed Stainless Steel Scoop,
12" X 6" X 2 3/4"
- 1101-20 Footed Polypropylene Scoop,
12" X 6" X 2 3/4"

SPECIFICATIONS

CAPACITY	
With Attachment Weights	2610g
Without Attachment Weights	610g
READABILITY	0.1g
BEAM CALIBRATIONS	
Center Beam	500g X 100g
Rear Beam	100g X 10g